






12.310 - Van Veen Grab – 2000 cm²

Manual



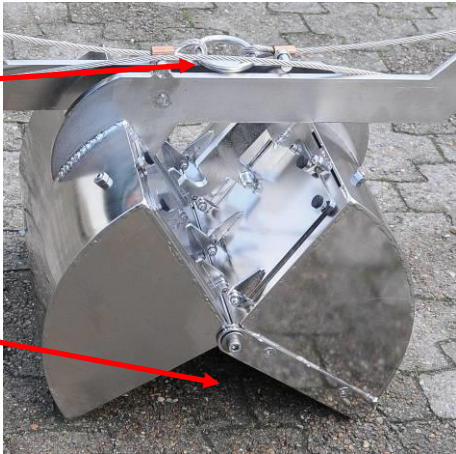


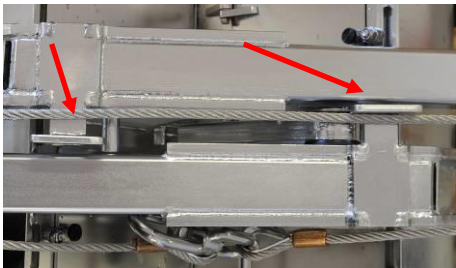
KC Denmark A/S






Research Equipment

Limnology • Oceanography • Hydrobiology

	<p align="center">Manual for Van Veen grab 2000 cm²</p>	<p align="center">Model no. 12.310</p>
	<p> Caution This Van Veen grab is very dangerous in unskilled hands and serious precautions must be taken to avoid accidents and injury to personnel.</p> <p>KC Denmark A/S is not, and cannot be held, responsible for any damage(s) made to equipment or to operators who ignore safety precautions or because of misuse or wrong operation.</p>	
	<p>Preparation:</p>	
<p>1</p>	<p> Caution</p> <p>Handle the grab with care to prevent injury to personnel. Fingers and hands may be stuck at the arms or by the jaws.</p>	
<p>2</p>	<p> Caution</p> <p>Inspect the thimbles and talurit ferrules for any damages or weakness. Replace if necessary. Otherwise, there may be a risk of losing the grab in the sea.</p>	

<p>3</p>	 <p>Caution</p> <p>Inspect the wire locks at both ends. It has to lock the wire to avoid the grab from hanging askew.</p>	
<p>4</p>	<p>Place the grab on a plain, stable surface. Force both arms down to horizontal position so the grab is open.</p>	
<p>5</p>	<p>For larger performance, you can add a maximum of 4 lead weights, 10 kg each (optional).</p> <p>Fasten the weights beneath the drainage holes. Use a M10 nut placed at the inner side of the grab.</p>	
<p>6</p>	<p>The locking pawl is shown in open (non-secured) position</p>	

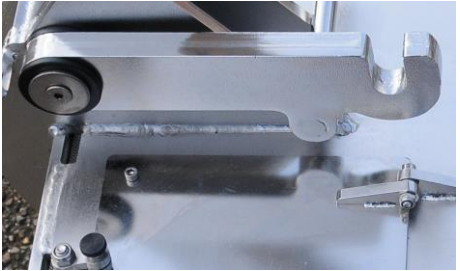7	Press the locking pawl upwards to lock the arms.	
8	From the upper side you can see if the pawl is locked properly. Now hold the steel wire straight to hold the pawl in its position.	
9	<p>The grab is ready and the wire for deploying the grab must be attached to the eye of the wire. The photo shows a smaller grab with thinner wire and accessories for lifting the grab.</p>  <p>Caution</p> <p>Keep the wires on the grab straight to avoid an unintended release of the locking mechanism.</p> <p>Recommended wire for deploying: Minimum 6-8 mm</p>	
10	 <p>Caution</p> <p>Align the wire at the inner side (towards the middle of the grab) of the two guiding rails; otherwise, the wire may get stuck at the arms upon release of the grab.</p>	

11	<p>Lower the grab to the sea bed at a reasonable speed. When the grab hits the sea bed, the wire will slacken and the locking pawl is released. Upon retrieving the grab, the shovels will close, and the sample will be taken.</p>	
Emptying the grab		
12	<p>The grab has 2 lids and turning the knob counter clockwise, the appropriate lid can be open and you can take out a small sample without emptying the whole sediment sample.</p> <p>Emptying the whole content of the grab is done by forcing the arms of the grab into horizontal position.</p>  <p>Caution</p> <p>Each time you handle the grab, take care to avoid any injury. See item #1</p>	
Maintenance		
13 A	<p>All parts of the Van Veen grab can be rinsed using salt water or fresh water. Regular cleaning with fresh water is recommended and all moveable parts must be moved individually to ensure all dirt has been removed.</p> <p>A. The locking pawl for the releaser and the hinges.</p>	
13 B	<p>B. The sieves in the drainage holes must also be free of sediment.</p>	
13 C	<p>C. The wheels at the end of the arms.</p>	

Technical information	
Materials:	
All metal parts:	AISI 316 stainless steel. Finish: Electro polish.
Wheels for wire:	Brass
Wire:	Stainless steel wire
Weights:	Lead, 10 kg each. Maximum mounting: 4 pcs.
Dimensions and weight:	
Overall dimensions, open, L x W x H:	245 x 55 x 63 cm
Overall dimensions, closed, L x W x H:	120 x 43 x 50 cm
Weight, exclusive lead weights:	95 kg net
Weight, inclusive 4 lead weights:	135 kg net
Sieves for the drainage holes:	Stainless steel, mesh size 1000 µm
Operational depth:	
Full oceanographic depth	>6000 m (no limits)

Rev. June 8, 2020 - lkj

KC Denmark A/S

Research Equipment
Limnology • Oceanography • Hydrobiology

E-mail: kc@kc-denmark.dk website: <http://www.kc-denmark.dk/>
 Holmbladsvej 19 - DK-8600 Silkeborg - Denmark - Tel. +45 86 82 83 47 - Fax +45 86 82 49 50
 Bank: Sydbank - S.W.I.F.T. SYBKDK22 IBAN DK5070460000104832 or
 Bank: Danske Bank - S.W.I.F.T. DABADKKK IBAN6730003462779860
 VAT no. DK 29 61 96 62